



qEVoriginal Rack


The **qEVoriginal Rack** is for the secure holding of qEVoriginal and qEV1* columns and sample collection tubes**. The rack is specifically designed with a tight fit to reduce movement during use. As a result, extreme care must be taken during assembly.

**qEV1 columns are compatible with the qEVoriginal Rack with the addition of the qEV1 Rack Accessory Pack.*

***1.5 mL sample collection tubes and 15 mL waste tubes are not provided. qEVoriginal and qEV1 columns are available from store.izon.com*

Care:



- ▶ Wash with warm water and soap.
- ▶ Do not use ethanol on this product.

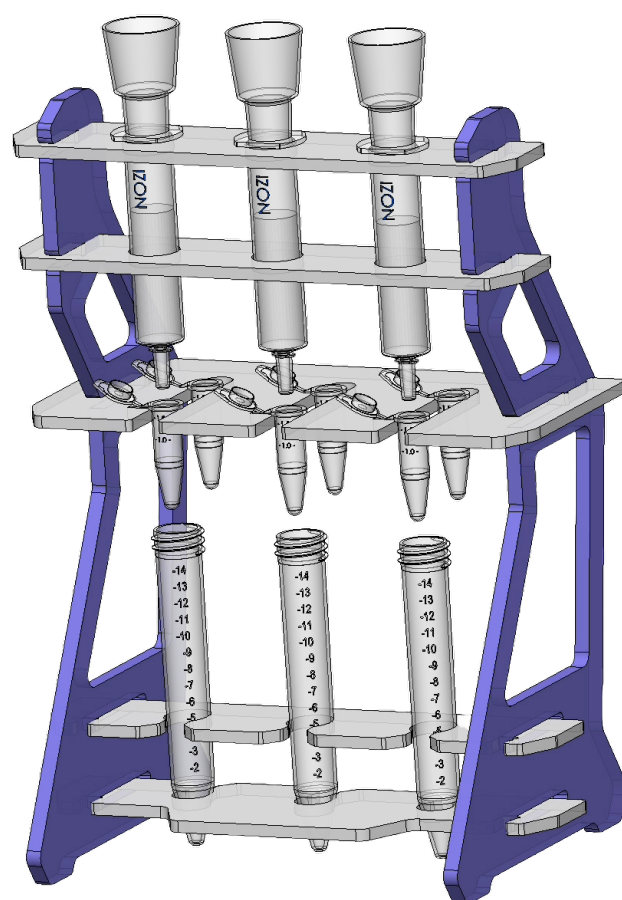

Avoid twisting joins or use of excess force during assembly, as this may result in the acrylic material cracking.


Once assembled, do not dismantle.

1

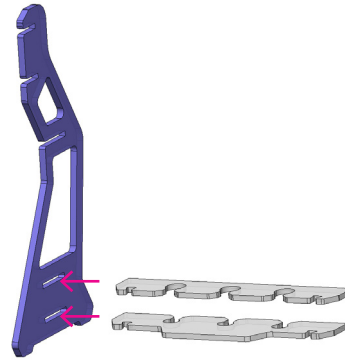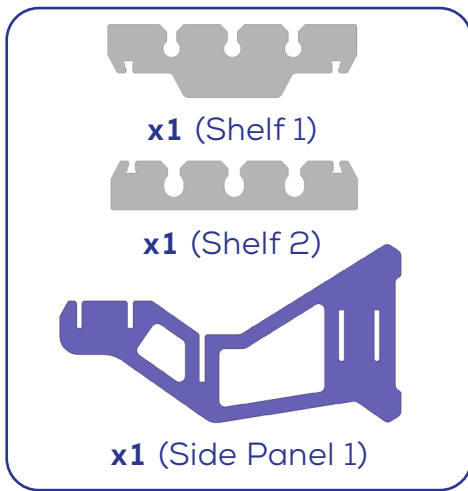



Begin by inserting one end of **Shelf 1** through the bottom slot of **Side Panel 1**, followed by **Shelf 2** in the slot above. Ensure both shelves click into place.

2

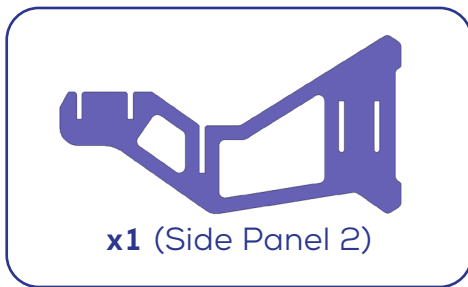


Insert the opposite ends of **Shelf 1** and **Shelf 2** into **Side Panel 2**. Once they have clicked into place, the rack should be freestanding.

3

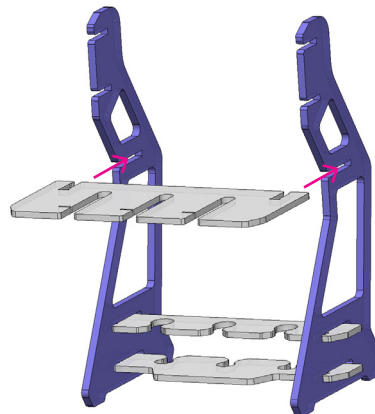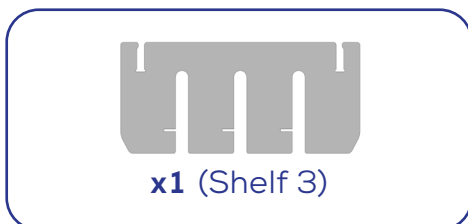

Position **Shelf 3** into the third slot in both **Side Panel 1** and **Side Panel 2**, ensuring it is secure.

4

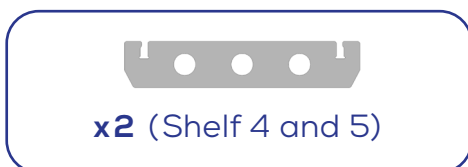

Insert **Shelf 4** and **Shelf 5** into the top two slots. Once all the shelves are secured, the rack is ready for use.