


qEVsingle Rack

The **qEVsingle Rack** is for the secure holding of qEVsingle columns and sample collection tubes*. The rack is specifically designed with a tight fit to reduce movement during use. As a result, extreme care must be taken during assembly.

**0.6 mL sample collection tubes and 1.5 mL waste tubes are not provided. qEVsingle columns are available from store.izon.com*

Care:



- ▶ Wash with warm water and soap.
- ▶ Do not use ethanol on this product.

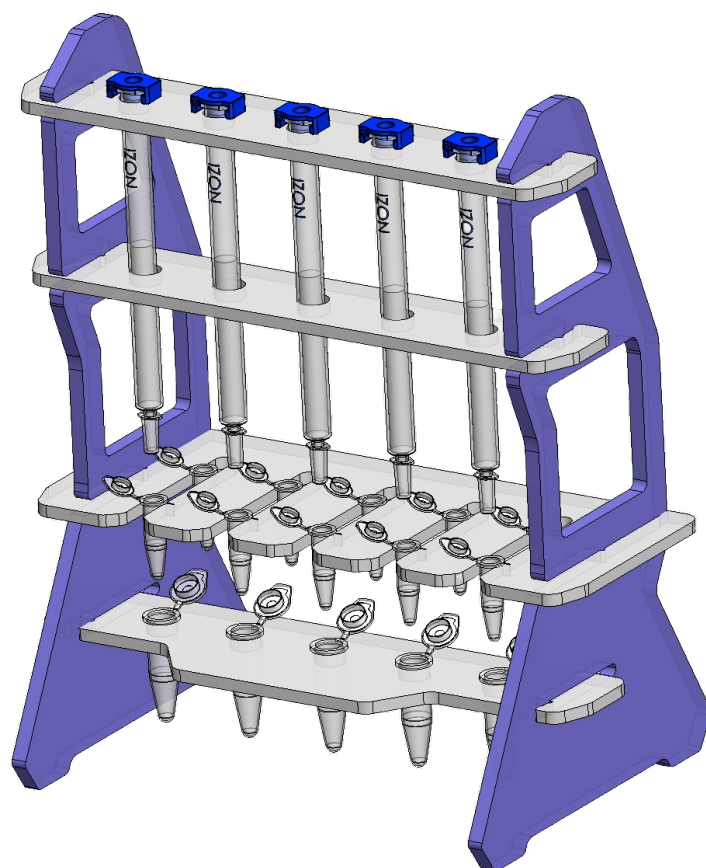

Avoid twisting joins or use of excess force during assembly, as this may result in the acrylic material cracking.



Once assembled, do not dismantle.

1

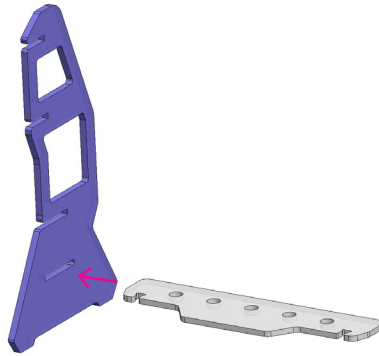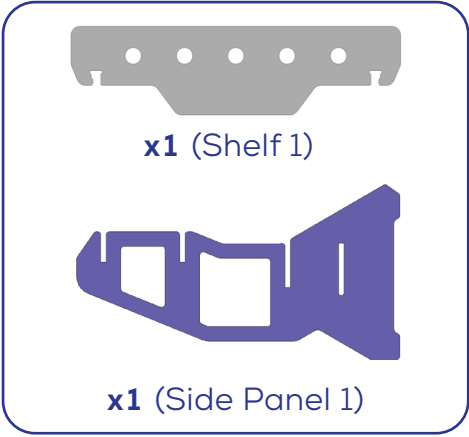


Begin by inserting one end of **Shelf 1** through the bottom slot of **Side Panel 1**. Ensure the shelf is clicked into place.

2

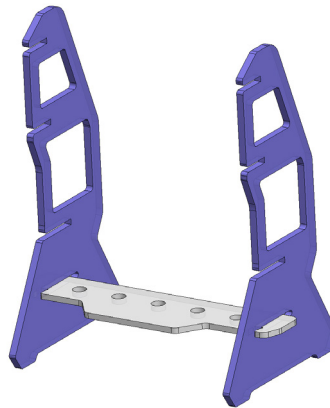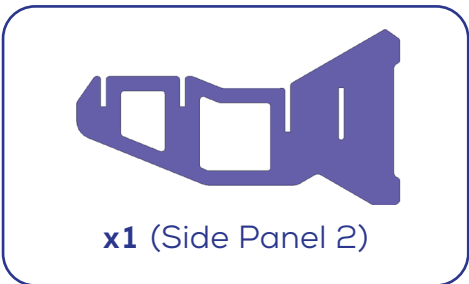



Insert the opposite end of **Shelf 1** into **Side Panel 2**. Once they have clicked into place, the rack should be freestanding.

3

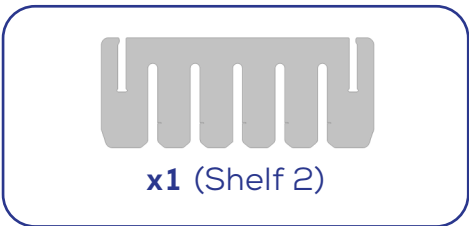

Position **Shelf 2** into the third slot in both **Side Panel 1** and **Side Panel 2**, ensuring it is secure.

4

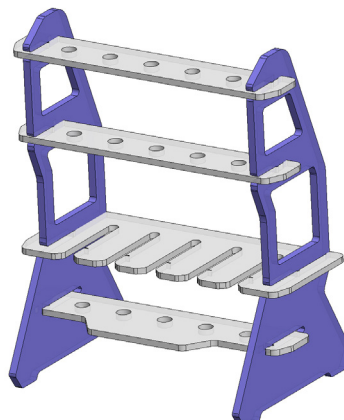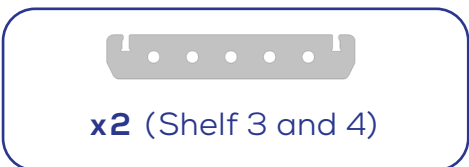

Insert **Shelf 3** and **Shelf 4** into the top two slots. Once all the shelves are secured, the rack is ready for use.